


# Norden Lab Compatibility Test

Some older Bioscreen C machines may not be compatible with Norden Lab Professional. Norden Logic developed a Compatibility Test program which is available from our web site free of charge. Follow the steps outlined below to install and run the test software successfully. The software can be found in the Bioscreen C download section at <http://www.nordenlogic.com>.


1. Download the installation ZIP file from Norden Logic.
2. Unzip the file into a folder of your choice.
3. Use the *Windows Explorer* or *Run...* to start the Compatibility Test program found in the folder where you unzipped the installation file (see example below).




4. Make sure all your Bioscreen C devices are connected and powered on.
5. Select *Search* to active automatic discovery of the attached Bioscreen C devices.

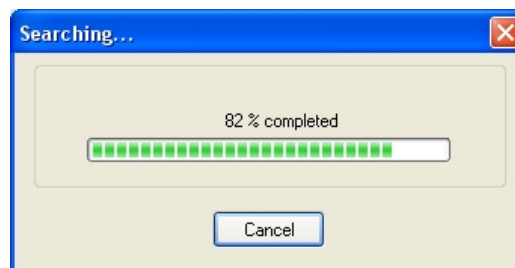


6. When the search is completed all found Bioscreen C devices are listed together with all RS-232 communications ports which do not have Bioscreen C devices attached to.

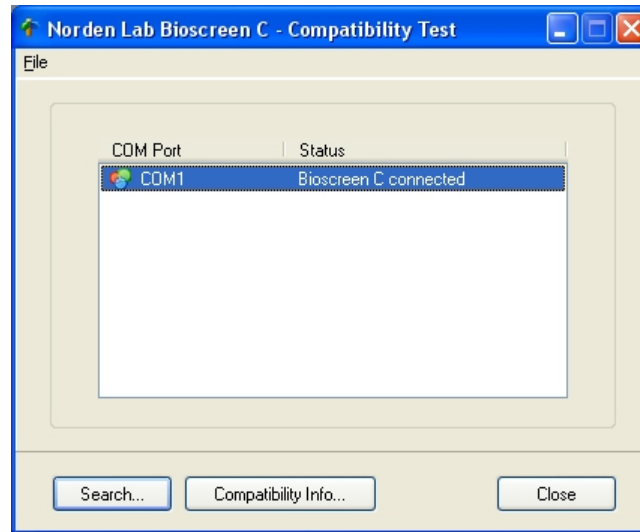


7. Select the Bioscreen C device you want to test. If a valid selection is selected *Compatibility Info* will be available. Click to view the information and to run the test.

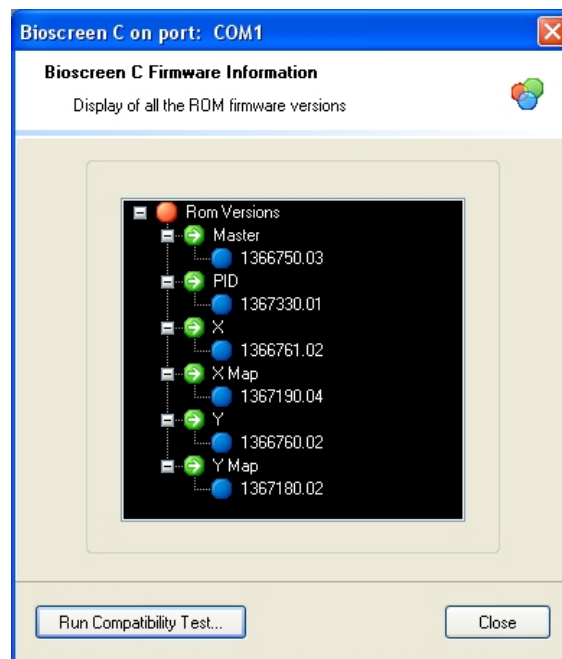

8. If you want to run the compatibility test click on *Run Compatibility Test* to bring up the test dialog box and click on *Start*.

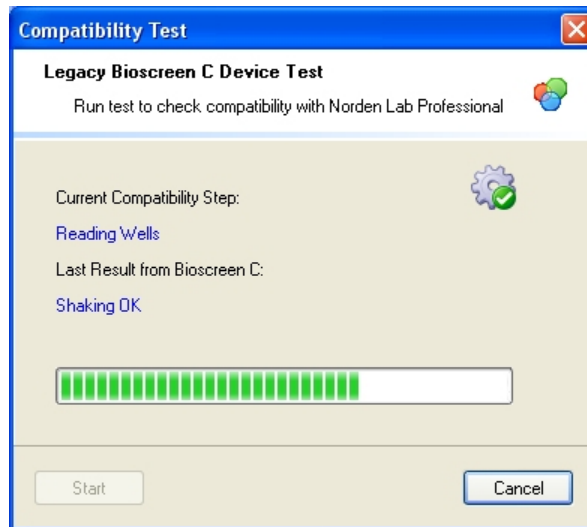

9. If the Bioscreen C device is 100% compatible with Norden Lab Professional a test result displaying a successful run will be shown. If an error was encountered during the run, please repeat the test. If the error persists, please contact your Bioscreen C distributor or Norden Logic for help.

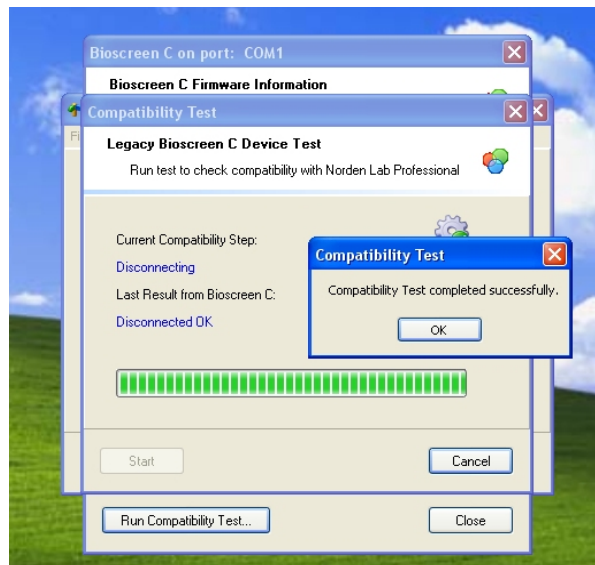

Revision 1.0A/1206