

March 2009

Windows Vista and Windows 7 have a security mode known as the User Account Control (UAC). This mode changes the role of the Administrator account. For more information regarding UAC please consult your Vista or Windows 7 online manual.

When installing Norden Lab Mastitis Studio General Edition for the first time you need to take the following into consideration in order to activate the 30 day evaluation period properly:

1. The user should be logged in as an Administrator
2. When starting the installation **right click** the installer and select "**Run as administrator**" and select "**Allow**" from the UAC control prompt.
3. After the installation completes the Administrator needs to run the application at least once as a true Administrator by **right clicking** the executable (from the menu "**Norden Lab Mastitis Studio General Edition**" or from folder (the file **MastitsStudioG.exe**) and select "**Run as administrator**" and select "Allow" from the UAC control prompt.

In case the software fails to activate correctly, please request a 30 day evaluation key from Finnzymes.

Note: Do not place the files into "C:\Program Files\...". Accept the installer default installation folder.

Note 2: For best Windows XP compatibility turn off the User Account Control (UAC) completely in the Accounts settings of Windows Vista or Windows 7. Please consult your administrator.